

REFRACTIVE INDEX LIQUID SERIES E

n(5893 Å) 25°C =1.5050

TYPICAL CHARACTERISTICS

COMPOSITION Triacetin and Ethyl Cinnamate
APPEARANCE Colorless to light yellow liquid
ODOR Fruity
COLOR STABILITY In sun: may slightly darken after 4 days; slightly more after 1 month; no significant change then after 10 years
INDEX CHANGE RATE BY EVAPORATION . Moderate: -0.00045 to -0.00074 expected: exposed surface area to volume ratio of 0.2 cm²/cc @ 25°C for 32 days
POUR POINT °C <7
BOILING POINT °C @ 760mm Hg >259
FLASH POINT °C COC >130
DENSITY g/cc @ 25°C 1.089
DENSITY TEMP. COEF. g/cc/°C -0.0010
COEF. OF THERM. EXP. cc/cc/°C 0.0009
VISCOSITY centistokes @ 25°C 12 (ca. 16 @ 15°C, 9 @ 35°C)
SURFACE TENSION dynes/cm @ 25°C .. 36
SOLUBLE: Acetone, Carbon Tetrachloride, Ethanol, Ethyl Ether, Freon TF, Methylene Chloride, Toluene, Xylene
PARTLY SOLUBLE: Heptane, Naphtha, Turpentine INSOLUBLE: Water
COMPATIBLE 3 week immersion @ 25°C: Mylar, Nylon, Polyester, Polyethylene, Polypropylene, Polyurethane, Phenolic, Teflon; Latex, Silicone, and Fluorosilicone Rubbers; Aluminum, Steel; (tests done on one example of each)
INCOMPATIBLE: Acrylic, Cellulose Acetate, Epoxy, Polycarbonate, Polystyrene, Polyvinyl Chloride; Neoprene; Tygon; may tarnish Copper and Brass
GEL FORMATION: may gel with age (very rare): 2 1/2 year shelf life
TOXICITY Low (request MSDS)

CAUCHY EQUATION: refractive index as a function of wavelength at 25°C

W = wavelength in angstroms (Å)

$$n(W) = 1.482607 + (583524.6)/W^2 + (6.736749E+12)/W^4$$

SOURCE OR SPECTRAL LINE	WAVELENGTH (angstroms)	REFRACTIVE INDEX 25°C	% TRANSMITTANCE 25°C		
			1mm	1cm	10cm
N laser	3370	1.586	75	6	0
i (Hg)	3650	1.564	99	90	36
h (Hg)	4047	1.5434	100	96	67
F' (Cd)	4800	1.5206	100	99	90
F (H)	4861	1.5194	100	99	91
e (Hg)	5461	1.5098	100	100	97
D (Na D1,D2 mean)	5893	1.5050	100	100	98
HeNe laser	6328	1.5014	100	100	98
C' (Cd)	6439	1.5006	100	100	98
C (H)	6563	1.4998	100	100	98
Ruby laser	6943	1.4976	100	100	99
GaAs laser	8400	1.4922	100	100	98
Nd:YAG laser	10648	1.488	100	98	82
Diode	13000	1.486	99	94	55
Diode	15500	1.485	99	88	27

$n_F - n_C$ = 0.0196
 Abbe v_D : $(n_D - 1)/(n_F - n_C)$ = 25.8
 Temp. coef: dn_D/dt 15-35°C = -0.000465

CARGILLE LABORATORIES INC.

55 COMMERCE RD. CEDAR GROVE N.J. 07009-1289 U.S.A.

phone: (973)239-6633 Fax: (973)239-6096